


Specification

model	WSC-2900 Rotator atto
speed	10 ~ 80rpm
angle	0°~ 90°
timer	1 ~ 9999min
power	AC100V 50/60Hz 35W
dimension	206(W) × 260(D) × 300mm(H)
weight	4.5kg

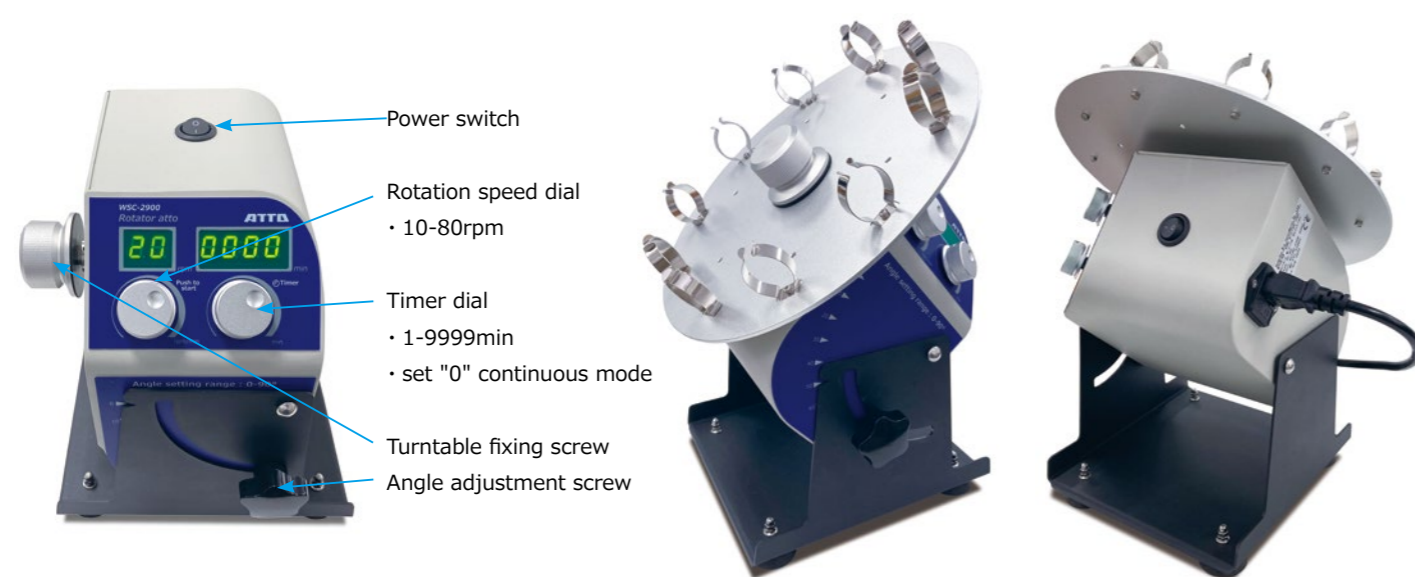

code	model	
4002900	WSC-2900 Rotator atto	1
4002940	Tunable A (for WSC-2900)	1
4002941	Tunable B (for WSC-2900)	1
4002942	Tunable C (for WSC-2900)	1
4002944	Tunable E (for WSC-2900)	1
4002949	Double platform connector (for WSC-2900)	1

Tabletop small volume rotary mixer

WSC-2900 Rotator atto

50ml conical tube
15ml conical tube
1.5ml eppendorf tube



Speed : 10 ~ 80rpm
Timer : 1 ~ 9999min
Angle : 0 ~ 90°

アトー株式会社



主要製品

- 発光・蛍光イメージングシステム
- 画像解析ソフトウェア
- 電気泳動装置
- 電気泳動関連試薬
- ウエスタンブロット試薬
- ペリスタポンプ
- 細胞培養・観察システム

- 東京本社 〒111-0041 東京都台東区元浅草3-2-2 ☎(03)5827-4861(代表) ☎(03)5827-6647
- 大阪支店 〒530-0044 大阪市北区東天満2-8-1 ☎(06)6136-1421(代表) ☎(06)6356-3625
若杉センタービル別館 5F
- 技術開発センター 〒110-0016 東京都台東区台東2-21-6 ☎(03)5818-7560(代表) ☎(03)5818-7563
◆メンテナンスサービスグループ ☎(03)5818-7567(代表) ☎(03)5818-7563

■URL <https://www.atto.co.jp/> お問い合わせ WEB会員登録の上お問い合わせフォームをご利用ください。



Rotator atto

Rotator atto is a rotary mixing apparatus. Mixing of 1.5ml / 15ml / 50ml tubes is possible.

Applications

- Extraction of samples such as DNA and proteins
- Culture mixing
- Immunoprecipitation


Function

- Angle : 0 - 90°
- speed : 10 - 80 rpm
- timer : 1 - 9999min
- timer[0] : contious operation



Angle setting

The angle of the turntable can be adjusted from the vertical state (0 °) to the horizontal state (90 °). By changing the angle, the solution in the tube can be agitated efficiently.

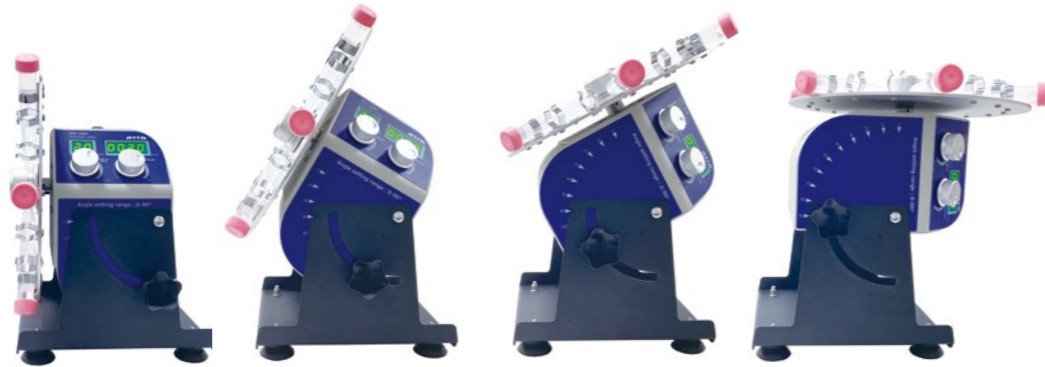

Stacking two turntables


Double platform connector for stacking two turntables(A/B/C/E). You can set various tubes together.



Option

Turntable A

Description 50ml tube × 10
Set the tube in a well-balanced manner.

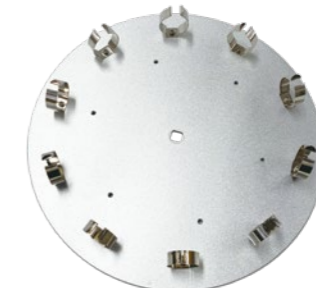

Turntabel A



Tube is sold separately

Turntable B

Description 15ml tube × 18
Set the tube in a well-balanced manner.




Turntabel B



Tube is sold separately

Turntable B

Description 1.5ml tube × 64
Set the tube in a well-balanced manner.

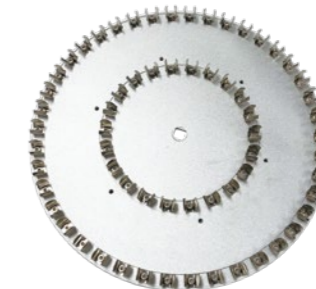


Turntabel C



Tube is sold separately

Turntable E

Description 50ml tube × 10
15ml tube × 10
1.5ml tube × 30
Set the tube in a well-balanced manner.

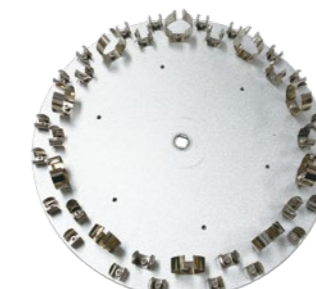

Turntabel E



Tube is sold separately

Double platform connector

Description Connector for stacking two turntables
for Turntable A/B/C/E



Double platform connector

